

Electronic Macro Nozzle Replacement Instructions



When to Replace:



If a nozzle is leaking or does not aspirate, a complete nozzle can be replaced.

Catalog No. 9060-5010

For electronic macro volume model 100–5000 μ L

Package includes:

- 2 nozzles
- Instructions

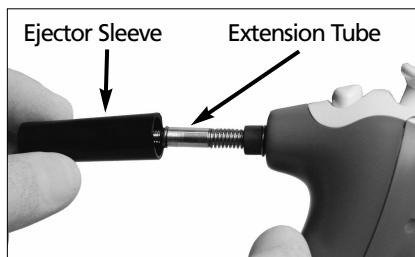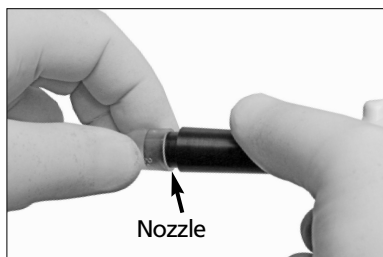

NOTE: Wear gloves when doing this procedure.

1. Place a macro pipette tip onto the nozzle and push on ejector sleeve until click is heard. Grip tip firmly near the nozzle, and turn it in a counterclockwise direction to unscrew and remove nozzle from pipette.
2. Slide off the ejector sleeve. Place a gloved finger lightly against the end of the air tube, and press and release plunger to expel any liquid that may be present in the end of the tube. Gently wipe the end of the air tube with a slightly dampened, lint-free tissue, then dry it off.
3. Replace the ejector sleeve. Note: The ejector sleeve is slightly narrower on the end that is closest to the pipette body.
4. Screw a new nozzle onto the extension tube. Tighten by hand, or use a macro pipette tip to ensure the nozzle is firmly attached to the extension tube.