

Please Read

When you first receive your ali-Q™, press the purple aliquoting button to wake the unit out of shipping mode.

For standard serological pipetting



Note: The motor will activate at seemingly random times when pressing the aspirate or dispense buttons. Don't worry. It's normal! For more information on ali-Q's intelligent measuring system, see Operating Manual at www.vistalab.com/support

For aliquoting



1. **Set aliquot volume** on dial
2. Aspirate adequate volume for total number of desired aliquots.



3. **Press & hold aliquot button** until the aliquot is dispensed. You will hear and feel the motor come on. Repeat for each aliquot. (For maximum precision, discard first aliquot.)

Note: if you don't hold the button long enough for the full aliquot volume, you will hear an "alert" sound and red light will flash.

**Description**

ali-Q pipet controller
ali-Q LS pipet controller

VistaLab Catalog #

2000-0005
2000-0003

Additional box contents:

spare 0.45 μ m membrane filter
power supply w/outlet adapters
wall mount
calibration certificate

9020-4005
9020-9005
9020-9008

**Register your ali-Q
online at
www.vistalab.com
for 1 year Warranty**

**Go to www.vistalab.com/support
for full Operating Manual**



Complete your optimal liquid handling system with	VistaLab Catalog #
Wobble-not™ serological pipets	
5 mL (200/cs)	4090-0005
10 mL (200/cs)	4090-0010
25 mL (200/cs)	4090-0025
50 mL (100/cs)	4090-0050



Think Differently. Feel Better.

2 Geneva Road Brewster, NY 10509 USA
(888) 652-6520; (914) 244-6226
customer.service@vistalab.com

©2017 VistaLab Technologies, Inc. All rights reserved.

L080-0052-001 RevB