¡IMPORTANTE! Lea este documento

Pipeteo serológico convencional

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LOB0-0052-001 RevC1

Register your ali-Q 2 online at www.vistalab.com for 1 year Warranty

Go to www.vistalab.com/support/ali-q-2 for full Operating Manual

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IMPORTANT! Please Read

Conventional serological pipetting

When you first receive your ali-Q™ 2, press the purple button to wake it out of shipping mode.

aspirate button
dispense button

Note:

1. The motor will activate at seemingly random times when pressing the aspirate or dispense buttons. Don't worry. It's normal!
2. ali-Q 2 will go into sleep mode after 3 minutes of non-use but will wake with any movement of the unit.

For more information on ali-Q's intelligent measuring system, see Operating Manual at www.vistalab.com/support

Aliquoting/repeat pipetting

speed dial

ali-Q 2 VS model only:
Set desired aliquoting speed using speed dial

aliquot dials

ali-Q 2 & ali-Q 2 VS ali-Q 2 LS

aliquot button

All ali-Q 2 models:

1. Set aliquot volume on aliquot dial.
2. Aspirate adequate volume for total number of desired aliquots.
3. Press and HOLD the purple aliquot button until the aliquot is dispensed. You will hear and feel the motor come on. Repeat press and hold for each aliquot. For maximum precision, discard first aliquot.

Note: if you don't hold the button long enough for the full aliquot volume, you will hear an "alert" sound and red light will flash.

Description

ali-Q 2 pipet controller	2100-0005
ali-Q 2 LS pipet controller	2100-0003
ali-Q 2 VS pipet controller	2100-1005

VistaLab Catalog

Additional box contents:

spare 0.22µm membrane filter	9020-4003
power supply w/outlet adapters	9020-9005
wall mount	9020-9008
calibration certificate	

Complete your optimal liquid handling system with: Wobble-not™ serological pipets VistaLab Catalog

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©2020 VistaLab Technologies, Inc. All rights reserved.
©2020 VistaLab Technologies, Inc. Tous droits réservés.