

The Ovation QS pipette is an adjustable volume, air displacement, 2 stroke piston pipette for aspirating and dispensing precise volumes.



Using your Ovation pipette





Ergonomic Handling

1. "Shake hands" position to pick up. Allow body to fill your palm.
2. Rotate adjustable hook to rest comfortably on your forefinger.
3. Keep hand posture relaxed.



Volume Setting

1. Press the center button  to cycle through 5 pre-set volumes, or press the up and down arrows  until desired volume appears.
2. "Lock in" volume by holding down plunger until three beeps are heard.



Installing the "Click Tip"

1. Insert nozzle into tip.
2. Lightly press down until a "click" is heard. This indicates the tip is securely sealed onto the nozzle.

Taking Care of Your Ovation Pipette

The Ovation pipette requires minimal routine maintenance. Always store it in its "standing" position when not in use. Clean outer surfaces as needed with a soft cloth dampened with warm water. To decontaminate outer surfaces, wipe with 70% ethanol or isopropanol solution or use a 10% bleach solution followed by water. Remove ejector sleeve and wipe outside of nozzle and inside of ejector sleeve with 70% alcohol to remove any accumulated deposits. **ONLY WIPE THE LCD SCREEN DISPLAY WITH WATER.**

- Performance specs guaranteed with VistaLab brand pipette tips only.
- 2 Year Limited Warranty.
- VistaLab Technologies offers comprehensive calibration and repair service. Annual service is recommended. See www.vistalab.com/services for more information or to place an order.

For more information on troubleshooting and maintenance visit www.vistalab.com/support.

For additional assistance with warranty repairs or other technical assistance, contact techservice@vistalab.com.



VistaLab Technologies, Inc.
2 Geneva Road
Brewster, NY 10509 USA
(888) 652-6520 or (914) 244-6226

Think Differently. Feel Better.