


VistaRak™ Autoclave Instructions

1. To prepare a VistaRak(s) for autoclaving, first remove all outer plastic wrapping.
2. Prop lid open using the locking tab.
3. Place VistaRak(s) in the autoclave at 121°C for 15 minutes, followed by a 15 minute drying cycle.
4. Close the cover before removing VistaRak(s) from autoclave to prevent any contamination.

Tips may be used after they have cooled to room temperature.

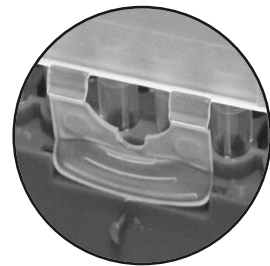


close up

Notes:

- Only non-filtered tips can be autoclaved.
- Ensure the VistaRak(s) have the proper bottom cover for autoclaving.

OK:

Flat, opaque bottom cover with recycle symbol

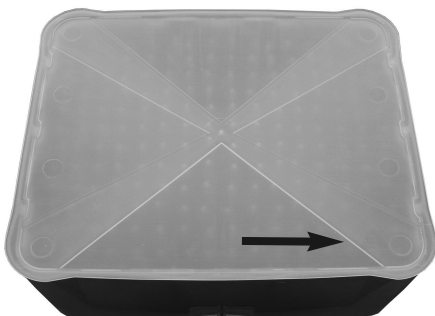

DO NOT AUTOCLAVE:

Clear older style bottom cover with ridges

